



How to return to the default settings


1. To restore the default settings of the camera we will go to the “**advanced settings**” of the software on the top right hand side of the screen.

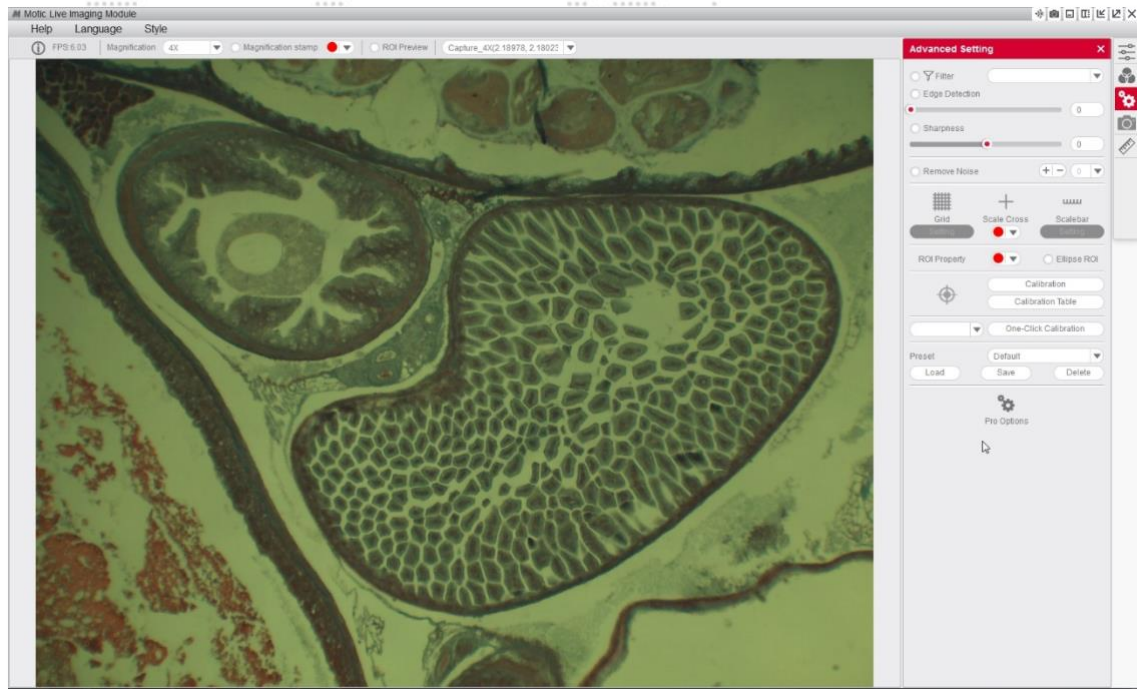


2. You will see a drop-down menu in the option “**Preadjust**” at the bottom of the screen. On this drop-down menu there is an option called “**Predetermined**”.

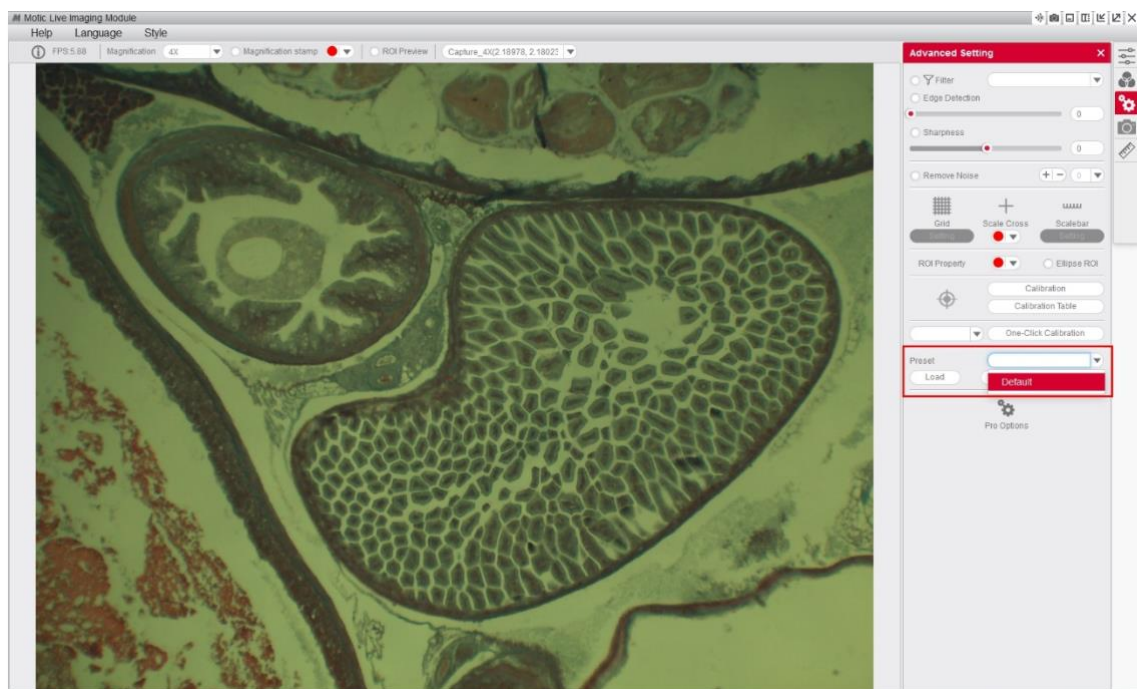

3. We will select the option “predetermined” and we will click on “Load”.

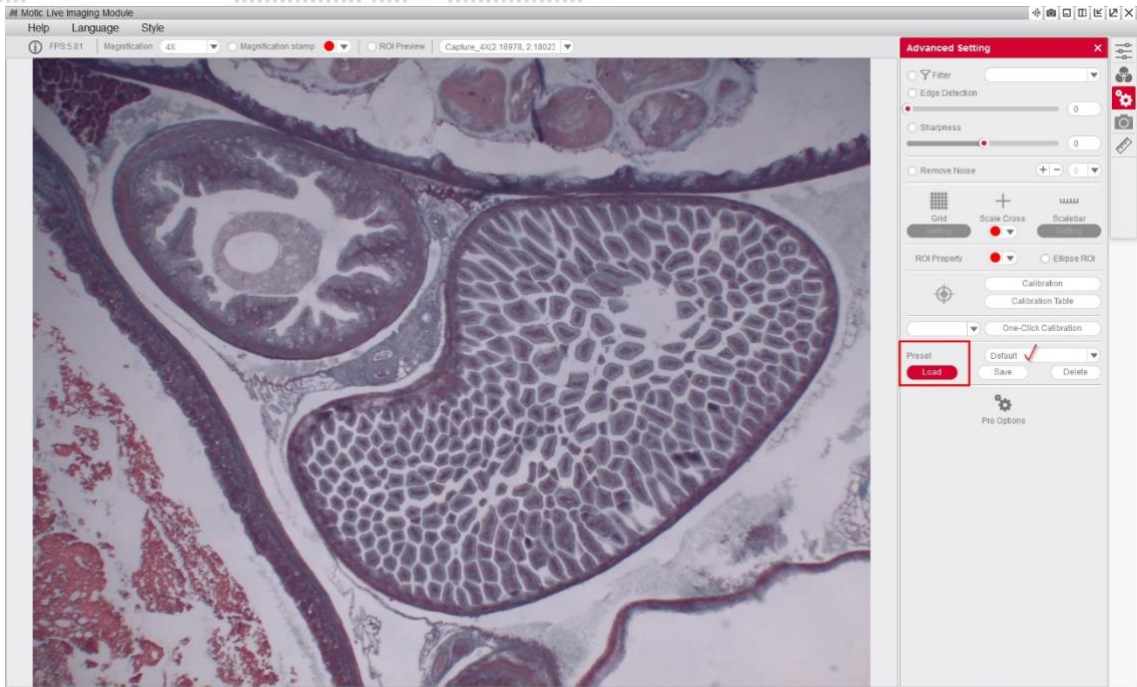

4. Now the software will have loaded the default settings of the camera.

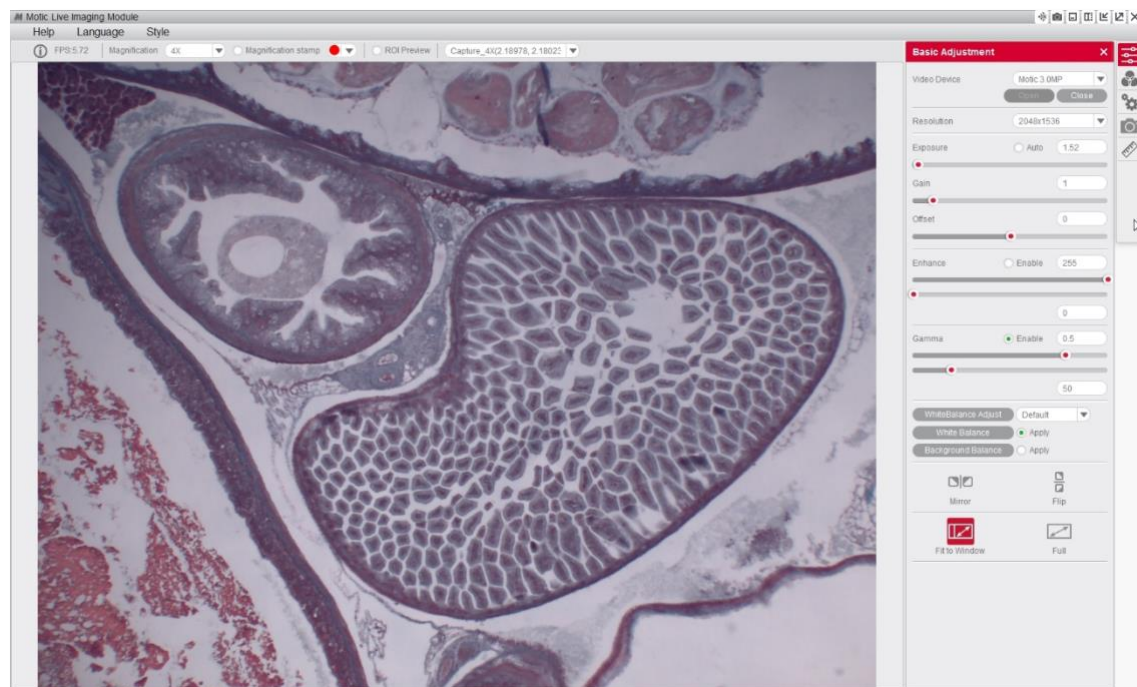

For more quick start guides to learn the different functions of our software, visit our web www.moticamseries.com

You can also contact our Digital Support Department by writing an e-mail to the following address: digital.support@moticeurope.com