
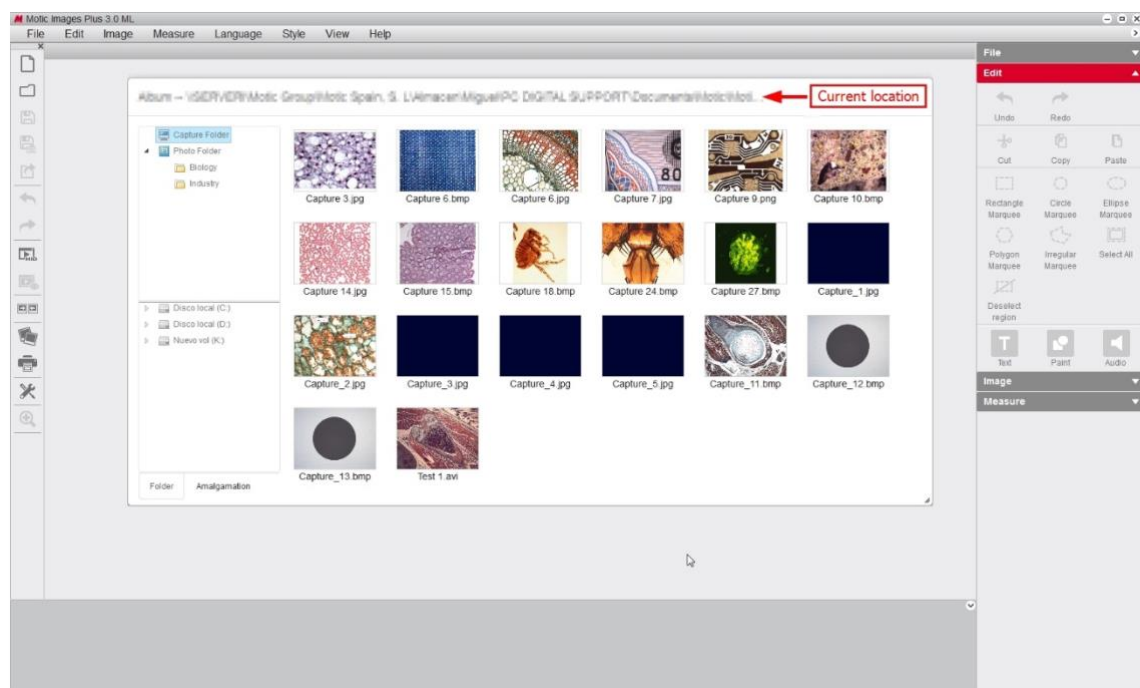


## How to relocate the Capture folder


All of the images that we capture with the software Motic Images Plus 3.0 are saved automatically in the “Capture folder” (default folder). This folder is created automatically when we install the software.

We can relocate the location of the captured images by choosing our preferred destination folder.

1. To do this, we need to go to the main screen of the software and select **“File-Album”**

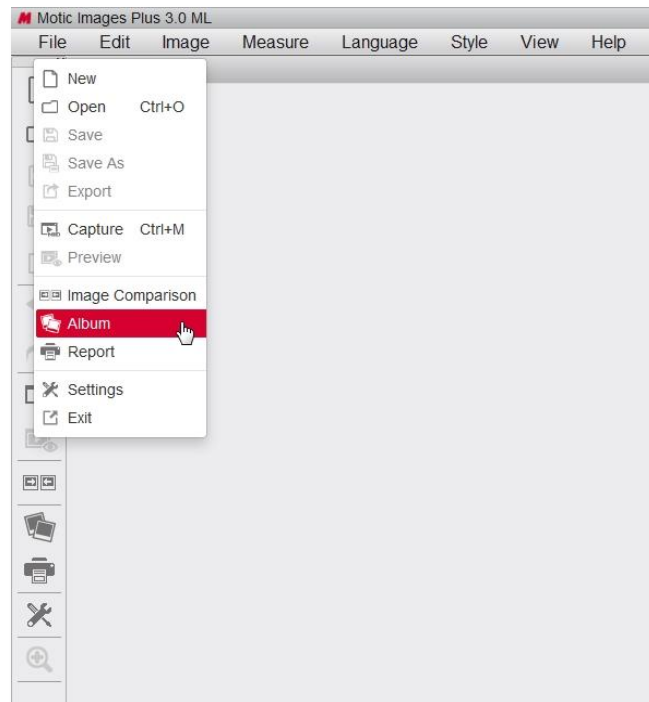


2. Right click with your mousebutton "Capture Folder". A menu will appear and here you will have to select "Set default capture folder".

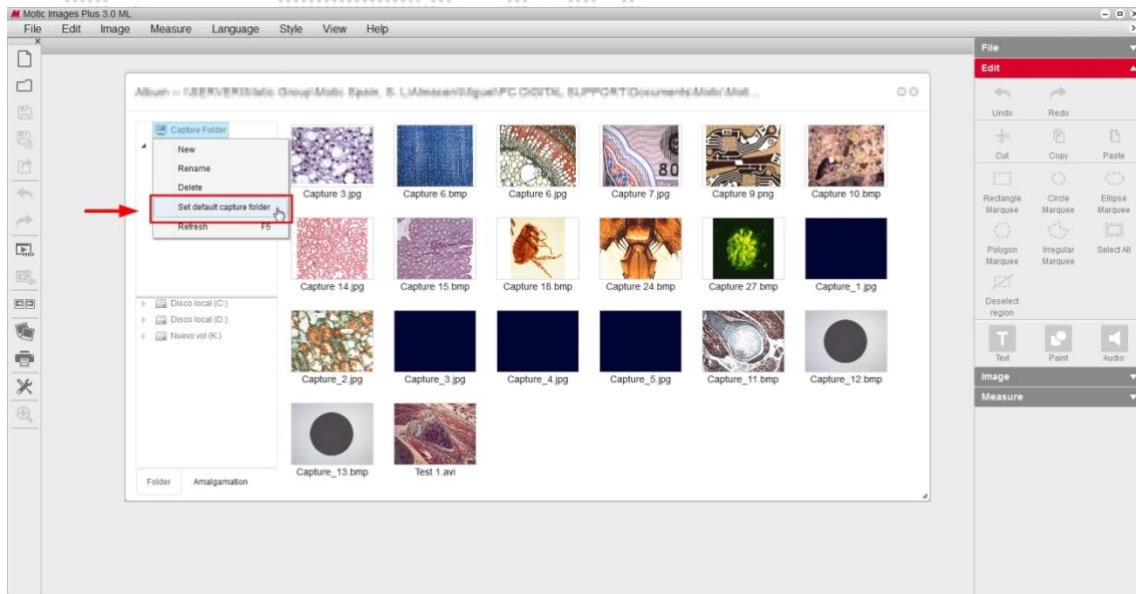

3. Here we will need to select the destination folder where we want the images to be saved automatically.

**NOTE:** We need to select a folder with write permissions in order for the images to access it.

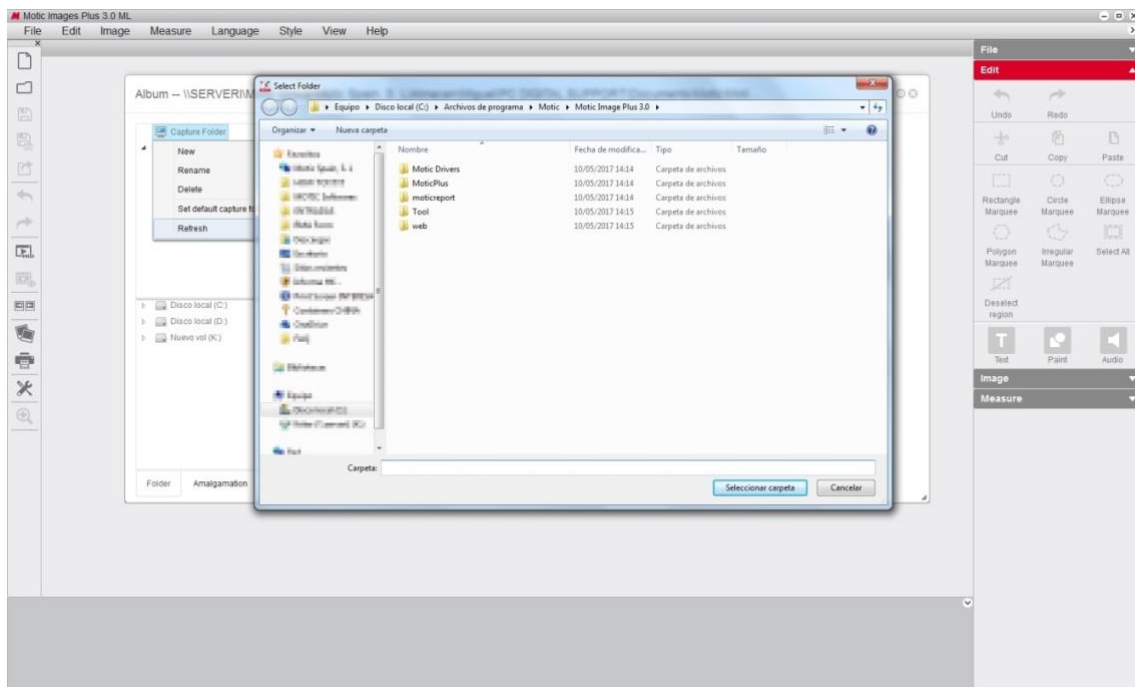

For more quick start guides to learn the different functions of our software, visit our web [www.moticamseries.com](http://www.moticamseries.com)

You can also contact our Digital Support Department by writing an e-mail to the following address: [digital.support@moticeurope.com](mailto:digital.support@moticeurope.com)