


How to choose the correct objective in the calibration table (live measurements)

***NOTE:** After calibrating the software (FAQ 2 How to calibrate with a dot) we will have to choose in the calibration table the objective that we have previously calibrated. For example: **4X-Motic**

1. Once we have calibrated and make the first measurement, we might see that the measurement is incorrect.



2. At the top of the software's tool bar we will see some objectives listed by default (which are not calibrated), as well as the new calibration that we ourselves have created. We need to select the same objective that we are using, and have previously calibrated, for example 4X-Motic. We will now see that the measurements are correct.

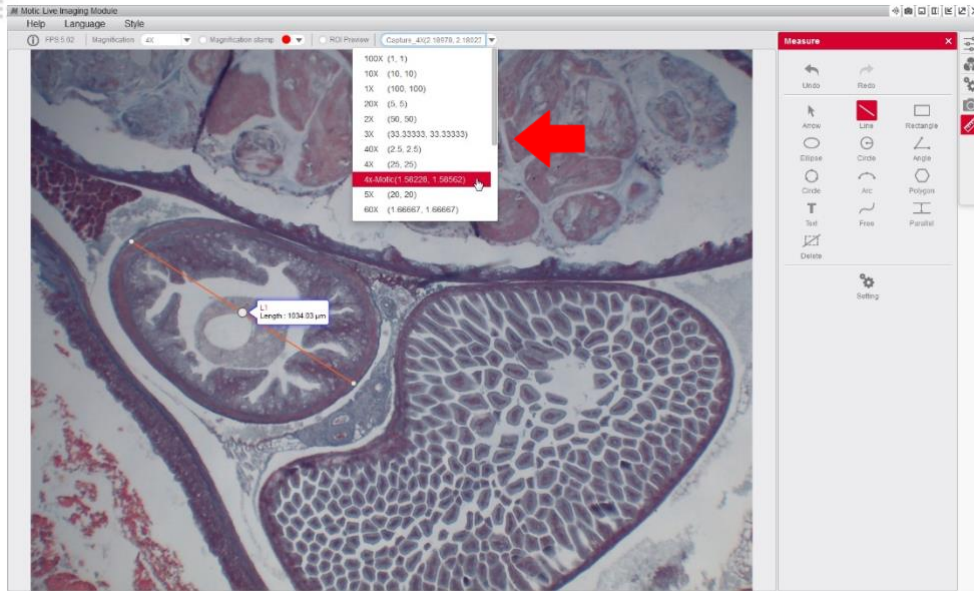

For more quick start guides to learn the different functions of our software, visit our web www.moticamseries.com

You can also contact our Digital Support Department by writing an e-mail to the following address: digital.support@moticeurope.com